

SCHEDULE 2

METHODS OF ANALYSIS

3a.

DETERMINATION OF NITRIC AND AMMONIACAL NITROGEN — ULSCH METHOD

1 SCOPE

1. This method is for the determination of nitric and ammoniacal nitrogen with reduction according to Ulsch.

2 FIELD OF APPLICATION

2. All nitrogenous fertilisers, including compound fertilisers, in which nitrogen is found exclusively in nitrate form, or in ammoniacal and nitrate form.

3 PRINCIPLE

3. Reduction of nitrates and nitrites to ammonia by means of metallic iron in an acid medium, and displacement of the ammonia thus formed by the addition of an excess of sodium hydroxide: distillation of the ammonia and determination of the yield of ammonia in a known volume of standard sulphuric acid solution. Titration of the excess sulphuric acid by means of a standard solution of sodium or potassium hydroxide.

4 REAGENTS

4

4.1 Hydrochloric acid solution, 50% (V/V): dilute an appropriate volume of hydrochloric acid ($d = 1.18 \text{ g/ml}$) with an equal volume of water.

4.2 Sulphuric acid, 0.1 N solution.

4.3 Sodium or potassium hydroxide, 0.1 N solution, carbonate free.

4.4 Sulphuric acid solution, approximately 30% H_2SO_4 (W/V), ammonia free.

4.5 Powdered iron reduced in hydrogen. (The prescribed quantity of iron must be able to reduce at least 0.05 g nitrate nitrogen).

4.6 Sodium hydroxide solution, 30 g per 100 ml, ammonia free.

4.7 Indicator solutions:

Mixed indicator:

(4.7.1) Solution A: dissolve 1 g methyl red in 37 ml 0.1 N sodium hydroxide solution and make up to 1 litre with water

Solution B: dissolve 1 g methylene blue in water and make up to 1 litre.

Mix 1 volume of solution A with 2 volumes of solution B.

This indicator is violet in acid solution, grey in neutral solution and green in alkaline solution; use 0.5 ml (10 drops).

Methyl red indicator solution:

Status: This is the original version (as it was originally made). This item of legislation is currently only available in its original format.

(4.7.2) dissolve 0.1 g methyl red in 50 ml 95% ethanol, make up to 100 ml with water and filter if necessary.

This indicator may be used (4-5 drops) instead of the preceding one.

4.8 Anti-bump granules of pumice stone, washed in hydrochloric acid and ignited.

4.9 Sodium nitrate.

5 APPARATUS

5. See Method 2.

6 PREPARATION OF SAMPLE

6. See Method 1.

7 PROCEDURE

7

Preparation of the solution

7.1 See Method 2.

Determination

7.2 Place in the receiving flask an exactly measured quantity of standard sulphuric acid (4.2) as indicated in the Table of Method 2 (variant (a)) and add the appropriate quantity of indicator solution (4.7.1 or 4.7.2).

The end of the extension tube of the condenser must be below, the surface of the standard acid in the receiving flask.

Using a pipette, transfer an aliquot part of the clear solution as indicated in the Table of Method 2 (variant (a)) and place it in the distilling flask of the apparatus. Add 350 ml water, 20 ml 30% sulphuric acid solution (4.4), stir, and add 5 g of reduced iron (4.5). Wash the neck of the flask with several ml water, and place in the neck of the flask a small, long stemmed funnel. Heat in a boiling water bath for an hour and then wash the stem of the funnel with a few ml water.

Taking care to avoid any loss of ammonia, add to the contents of the distilling flask 50 ml concentrated sodium hydroxide solution (4.6), or in the cases where 20 ml hydrochloric acid (4.1) has been used to dissolve the sample, add 60 ml of concentrated sodium hydroxide solution (4.6). Assemble the distillation apparatus. Distil the ammonia according to the procedure given in Method 2. Titrate the excess acid with the standard solution of sodium or potassium hydroxide (4.3).

Blank test

7.3 Carry out a blank test (omitting only the sample) under the same conditions and allow for this in the calculation of the final result.

Control test

7.4 Before analysis, check that apparatus is working properly and that the correct application of the method is used by using an aliquot part of a freshly prepared solution of sodium nitrate (4.9) containing 0.045 g to 0.050 g of nitrogen.

Status: This is the original version (as it was originally made). This item of legislation is currently only available in its original format.

8 EXPRESSION OF THE RESULTS

8. Express the results of analysis as a percentage of nitric nitrogen or combined ammoniacal and nitric nitrogen contained in the fertiliser as received for analysis.