
Status: EU Directives are being published on this site to aid cross referencing from UK legislation. After IP completion day (31 December 2020 11pm) no further amendments will be applied to this version.

[^{F1}ANNEX I

TEST SCHEME FOR DIAGNOSIS, DETECTION AND IDENTIFICATION OF THE RING ROT BACTERIUM, *CLAVIBACTER MICHIGANENSIS* (Smith) Davis *et al.* ssp. *SEPEDONICUS* (Spieckermann et Kotthoff) Davis *et al.* SCOPE OF THE TEST SCHEME

Textual Amendments

- F1** Substituted by [Commission Directive 2006/56/EC of 12 June 2006 amending the Annexes to Council Directive 93/85/EEC on the control of potato ring rot.](#)

Status: EU Directives are being published on this site to aid cross referencing from UK legislation. After IP completion day (31 December 2020 11pm) no further amendments will be applied to this version.

Appendix 3

Buffers for test procedures

1. Buffers for extraction procedure

1.1. Extraction buffer (50 mM phosphate buffer, pH 7,0)

This buffer is used for extraction of the bacterium from plant tissues by homogenisation or shaking.

Na ₂ HPO ₄ (anhydrous)	4,26 g
KH ₂ PO ₄	2,72 g
Distilled water	1.00 L

Dissolve ingredients, check pH and sterilise by autoclaving at 121 °C for 15 min.

Additional components may be useful as follows:

	Purpose	Quantity (per L)
Lubrol flakes	Deflocculant ^a	0,5 g
DC silicone antifoam	Anti-foam agent ^a	1,0 ml
Tetrasodium pyrophosphate	Anti-oxidant	1,0 g
Polyvinylpyrrolidone-40 000 (PVP-40)	Binding of PCR inhibitors	50 g

^a For use with homogenisation extraction method.]
